

Technical Data Sheet

1. Product


103.3C - pH Indicator Strips, 0 - 6.0.

2. Instructions for use

- Immerse an indicator strip in the test solution for 2 seconds.
- Shake off excess solution.
- Wait 10-20 seconds.
- Compare the colour produced on the strip against the colour chart before the indicator pads dry out.

3. Measuring range

pH 0.0 - 0.5 - 1.0 - 1.5 - 2.0 - 2.5 - 3.0 - 3.5 - 4.0 - 4.5 - 5.0 - 5.5 - 6.0

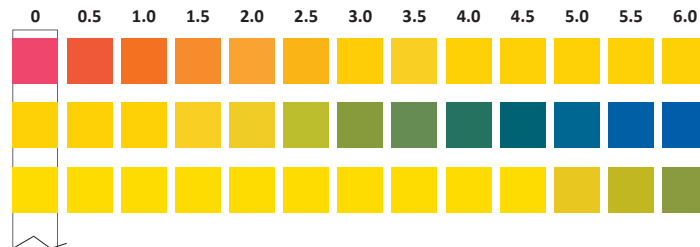

4. Application

pH indicator strip for testing of pH in solutions between pH 0.0 and pH 6.0.

5. Shelf Life

3 years from manufacturing date.

6. Contents

100 strips.

7. Weights & Dimensions

pH strip: 5m x 85mm

Plastic pack: 65mm(w) x 88mm(h) x 15mm(d)

8. Notes

Keep any unused strips in the original container. Do not store where they are liable to be exposed to acid or alkaline gases or vapours. Store in a cool dry place.